

**TECHNICAL SHEET**

<b>B600</b>	<b>MRS BROTH MODIFIED (LACTOBACILLI HETEROFERM SCREEN BROTH)</b>				
<b>Formula</b>					
<b>Ingredients :</b>		<b>gms/lit.</b>			
Dextrose		20.00			
Proteose peptone		10.00			
Yeast extract		5.00			
Sodium acetate		5.00			
2-Phenylethyl alcohol		3.00			
Ammonium citrate		2.00			
Dipotassium phosphate		2.00			
Magnesium sulphate		0.10			
Manganese sulphate		0.05			
Bromo cresol green		0.04			
Cycloheximide		0.004			
Final pH (at 25°C) :		4.3 ± 0.2			
<b>Directions :</b>					
Suspend 47.2 grams in 1000 ml distilled water containing 1 ml polysorbate 80. Mix thoroughly and dispense in tubes containing inverted Durhams tubes. Sterilize by autoclaving at 15 lbs pressure (121°C) for 15 minutes. If necessary, adjust the pH with glacial acetic acid after sterilization.					
Warning: Cycloheximide is very toxic. Avoid skin contact or aerosol formation and inhalation.					
<b>Principle :</b>					
Proteose peptone and dextrose supply nitrogen, carbon and other elements essential for the growth of Lactobacilli. Polysorbate 80 a mixture of oleic esters, supplies fatty acids required by Lactobacilli. Ammonium citrate, Sodium acetate, 2-phenylethyl alcohol and Cycloheximide inhibit gram – negative organisms, molds and certain gram – positive bacteria. Certain yeasts are suppressed because of presence of Cycloheximide.					
<b>QC Tests – (I) Dehydrated Medium</b>					
Colour :		Light yellow to bluish grey			
Appearance :		Homogeneous Free Flowing powder			
<b>(II) Rehydrated medium</b>					
pH (post autoclaving/heating) :		4.3 ± 0.2			
Colour (post autoclaving/heating) :		Green			
Clarity (post autoclaving/heating) :		Clear to slightly opalescent			
<b>(III) Q.C. Test Microbiological</b>					
Cultural characteristics observed after upto 3 days at 35-37°C with 5 – 10% CO <sub>2</sub> .					
MICROORGANISM (ATCC )		GROWTH			
Lactobacillus plantarum (8014 )		Luxuriant			
Lactobacillus fermentum (9338 )		Luxuriant			
Lactobacillus acidophilus (4356 )		Luxuriant			
Lactobacillus casei (9595)		Luxuriant			
<b>Precautions :</b>		1. For Laboratory Use.			
		2. Follow proper, established laboratory procedures in handling and disposing of infectious materials.			
<b>Limitations :</b>		1. Since the nutritional requirements of organisms vary, some strains may be encountered that fail to grow or grow poorly on this medium.			
<b>Use :</b>		For the isolation and cultivation of Lactobacillus species from salad dressings.			
<b>Storage :</b>		Dehydrated medium and prepared medium – Between 2 to 8°C.			
<b>Packing :</b>		500 gm. bottle			
<b>Product profile:</b>		Reconstitution	Quantity on Preparation (500g)	pH (25°C)	Supplement
<b>B600</b>	47.2g/l		10.593L	4.3 ± 0.2	Polysorbate 80
					121°C / 15 minutes

Refer disclaimer Overleaf

**BIOMARK Laboratories-INDIA**

[www.biomarklabs.com](http://www.biomarklabs.com)

**TECHNICAL SHEET**

User must ensure suitability of the product(s) in their application prior to use. Products conform solely to the information contained in this and other related BIOMARKLABORATORIES publications.

The information contained in this publication is based on our in-house studies and market performance and is to the best of our knowledge true and accurate. BIOMARK LABORATORIES reserves the right to make changes to specifications and information related to the products at any time. Products are not intended for human or animal or therapeutic use but for laboratory, diagnostic, research or further manufacturing use only, unless otherwise specified. Statements contained herein should not be considered as a warranty of any kind, expressed or implied, and no liability is accepted for infringement of any patents.

Page 02 of 02