

**BF129**

**ANTHERACIS SELECTIVE SUPPLEMENT**

**FORMULA**

**Ingredients :**

(Per vial, sufficient for 1000 ml medium)

Lysozyme	300000 Unit
Polymyxin B sulphate	30000 Unit

**Directions :**

Rehydrate the contents of 1 vial aseptically with 10 ml sterile distilled water. Mix well and add aseptically to sterile molten, cooled to (45-50°C) PLET Agar Base (B1242) / PLET Agar Base, Modified (B1243). Mix well and dispense as desired.

**Precautions :**

1. For Laboratory Use.
2. Follow proper, established laboratory procedures in handling and disposing of infectious materials.

**Use :** Anthracis Selective Supplement is recommended for the selective isolation of *Bacillus anthracis*

**Storage:** Between 2 to 8°C.

**Packing :** 5 vial